Digital Refrigerator Thermometer

TMX - Series

Index

	1
	1
TMX - Series	1
	1
Index	2
Overview	3
Safety Instructions	4
Please wear nitrile rubber gloves in case of a leaky bottle.	4
Unpacking	5
Delivery contents	5
.Implementing	6
Standard usage	6
Display	7
Buttons	7
Handling	8
What to do, if...	10
Repair and Service	10
Calibration	11
Warranty	11
Cleaning and Maintenance	12
Disposal	12
Technical Data	13
Approvals	14

Overview

The TMX thermometer series is for measurement of ambient temperature, but also to measure in cooling and freezing areas designed.

The following types are available:

TMX 310:

It has an internal temperature sensor and an external temperature probe which is in a bottle with glycol.

TMX 320:

It has an internal temperature sensor and an external temperature probe.

TMX 410:

It has two external temperature probes, each in a bottle of glycol.

TMX 420:

It has two external temperature probes.

Safety Instructions

Tightness:

The thermometer TMX 310 and TMX 410 are delivered with glass bottles of glycol solution.

Please check the bottles of possible existing cracks.

**ATTENTION HAND SAFETY,
CUT RISK!**

Please mind eventual leak of the glycol solution.

Please wear nitrile rubber gloves in case of a leaky bottle.

Unpacking

Please take all parts out of the box.

Carefully handle the sensor cables so that you do not damage them.

Check the glass bottles for possible damage and for leaking glycol during and after unpacking.

Delivery contents

The package includes the following parts:

- Thermometer
- Battery (Type AAA)
- Wall bracket
- Velcro® tape
- User Manual
- Certificate of Calibration

Implementing

1. Unpack all the delivered parts.
2. Check if the delivery is complete.
3. Put the battery in the device.
4. Choose the temperature unit between °C and °F.
5. Control the limit settings in the thermometer.
6. Activate the alarm function.
7. Put the thermometer on its final destination.
8. Place the probes on the right location. Make sure that the cable needs to be pinched in the door seal.

Standard usage

The Refrigerator Thermometer is for use in ambient conditions.

It is designed to control the temperature in cooling and freezing areas in laboratories and drug stores.

Display

The Thermometer shows current and Max/Min value at the same time.

Buttons

Alarm-Confirmation Button

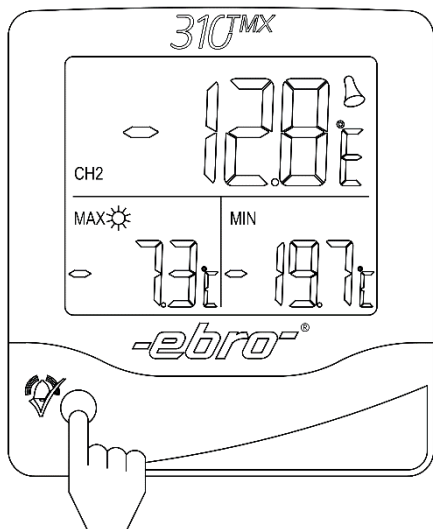
The Alarm-Confirmation button is placed on the front side of the device.

Following buttons are located on the rear side of the device:

Channel Switch

Alarm button

Button to choose the temperature unit

Handling

Alarm Function

The thermometer has an optical and acoustical alarm.

This function can be activated or even deactivated by pushing the alarm button.

Additionally to the standard alarm function, the acoustic alarm can be confirmed by pushing the Alarm Confirmation button while an acoustic alarm is active.

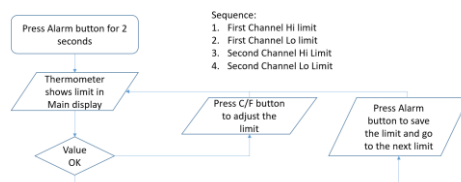
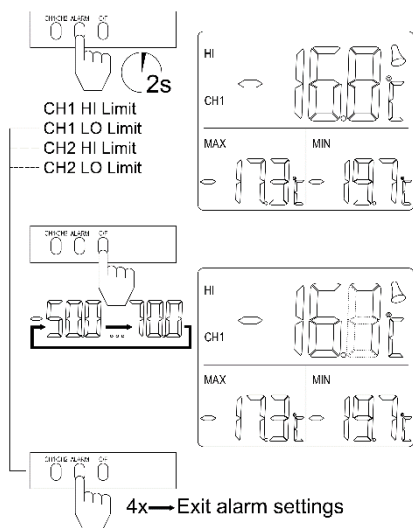
The optical alarm stays until the measured value is back between the limits.

Switch Alarm function on / off

Push the Alarm button for 1 second to turn the alarm function on or off.

Set Limits

Push the alarm button for 3 seconds to start limit settings. Follow the procedure.

Alarm Confirmation

Push the Alarm Confirmation button while the acoustic alarm is active to quit the alarm.

Clear Max/Min Memory

Push the Alarm Confirmation button for at least 3 seconds to clear the Max/Min memory.

Choose temperature unit

Push the temperature unit button to switch between °C and °F.

Insert battery

1. Fold out the stand.
2. Open the battery compartment by unlocking and moving the battery cover.
3. Now unpack the battery and place it in the battery holder. Pay attention to the polarity of the battery.
4. Close the battery compartment.

What to do, if...*No Display*

Check if the battery was inserted correctly.

Please replace the battery.

Display shows „Lo.“

The measured value is outside of the measurement range.

Display shows „Hi.“

The measured value is outside of the measurement range.

Display shows “err”

Probe or cable is broken.

No Alarm Signal

Alarm function wasn't activated.

Repair and Service

If you have any complaints about the TMX, please contact our customer service.

You will find additional information at:
www.ebro.com

Calibration

Each thermometer should be maintained and calibrated from time to time.

We don't offer any calibration service for thermometers from the series TMX.

We recommend to buy a new TMX thermometer instead of calibration.

The TMX is delivered with a Certificate of Calibration which is 2 years valid from the first calibration during the production.

Warranty

We grant a 2 year warranty on the function of the device.

Cleaning and Maintenance

Use a damp cloth to clean the TMX thermomixer.

Do not use a solvent, such as acetone, because this damages the plastic.

Disposal

If the device becomes no longer fit for purpose, it must be disposed of in a suitable, environmentally-friendly manner.

Technical Data

Operating Temperature	-20 °C ... +50 °C
Storage Temperature	-20 °C ... +60 °C
Temperature Units	°C, °F
Accuracy	±0.5 °C (-20 °C ... +40 °C) ±1.0 °C remaining area
Resolution	0.1 °C
Response time	30 seconds
Housing Material	ABS
Protection rate	IP 20
Dimensions	
(L x W x H)	100 x 110 x 23 mm
Cable length	3 m
Weight	app. 120 g
Battery	AAA (Alkaline)
Battery lifetime	app. 12 months, without alarms

TMX 310

Measuring range	
Internal Sensor	-20 °C ... +50 °C
External probe	-50 °C ... +70 °C
Probe type	glycol bottle

TMX 320

Measuring range	
Internal Sensor	-20 °C ... +50 °C
External probe	-50 °C ... +70 °C
Probe type	metal capsule 4 mm diameter, 20 mm length

TMX 410

Measuring range	-50 °C ... +70 °C
Probe type	glycol bottle
Amount of probes	2

TMX 420

Measuring range	-50 °C ... +70 °C
Probe type	metal capsule 4 mm diameter, 20 mm length
Amount of probes	2

Approvals

Operation is subject to the following two conditions:

- (1) This device may not cause harmful interference, and
- (2) This device must accept any interference received including interference that may cause undesired operation.

Changes or modifications of the device could void the user's authority to operate the equipment.

Users manual or instruction manual may downloaded at www.ebro.com over the internet.

Service-Adresse / Service Address / Adress du Service

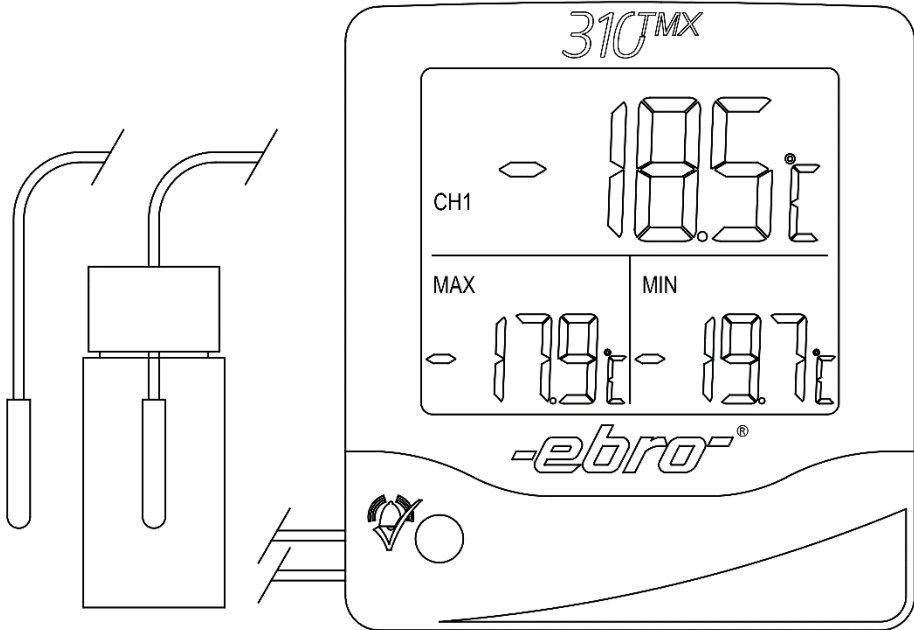
Xylem Analytics Germany Sales GmbH & Co. KG
ebro
Dr.-Karl-Slevogt-Str. 1
82362 Weilheim
Germany

Phone: +49.(0)841.954.78.0
Fax: +49.(0)841.954.78.80
Internet: www.ebro.com
E-Mail: ebro@xyleminc.com

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Xylem Analytics Germany GmbH